



## UNPACKING INSTRUCTIONS


# Bead Ruptor 96+

Version: A  
Document Number: 03-300

Date: August 2024

# Unpacking Instructions


**NOTE:** There is a ShockWatch indicator on the packaging box, which is used to check for improper handling of the shipment. Upon delivery, check if the indicator is red. In this case, follow the instructions on the ShockWatch label.

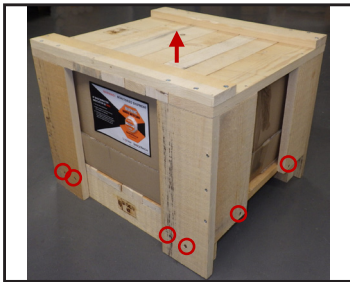

1. Unscrew 12 screws at the bottom of wooden crate. 6 screws are marked on above picture with circles, other 6 screws are located at the same positions on the other side of the crate. Then remove the upper part of the wooden crate by lifting it.



2. Remove the upper part of cardboard box by lifting it.

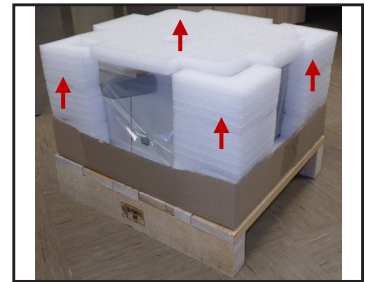


3. Remove the packaging materials.



4. The weight of the machine is 45.5 kg (100 lbs). To prevent possible injuries, at least two people should lift and carry the machine by holding it at the bottom from opposite sides. Take the machine out of the box and put it on the table.



5. Remove the plastic bag from the machine.

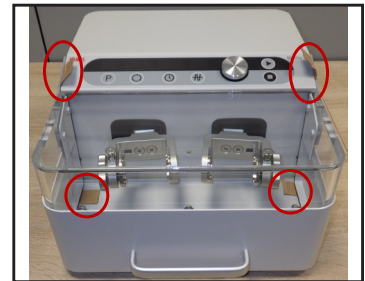

6. Open the lid and remove the cardboard pieces used for protection of the lid.



7. Close the lid and connect the power cord.



8. Check the operation of the machine.



The information provided in this document is for reference purposes only and may not be all-inclusive. Revvity, Inc., its subsidiaries, and/or affiliates (collectively, "Revvity") do not assume liability for the accuracy or completeness of the information contained herein. Users should exercise caution when handling materials as they may present unknown hazards. Revvity shall not be liable for any damages or losses resulting from handling or contact with the product, as Revvity cannot control actual methods, volumes, or conditions of use. Users are responsible for ensuring the product's suitability for their specific application. REVVITY EXPRESSLY DISCLAIMS ALL WARRANTIES, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, REGARDLESS OF WHETHER ORAL OR WRITTEN, EXPRESS OR IMPLIED, ALLEGEDLY ARISING FROM ANY USAGE OF ANY TRADE OR ANY COURSE OF DEALING IN CONNECTION WITH THE USE OF INFORMATION CONTAINED HEREIN OR THE PRODUCT ITSELF.

[www.revvity.com](http://www.revvity.com)

revvity

For research use only. Not for use in diagnostic procedures.

Revvity, Inc.  
940 Winter Street  
Waltham, MA 02451  
USA  
[www.revvity.com](http://www.revvity.com)

For a complete listing of our global offices, visit [www.revvity.com](http://www.revvity.com)  
Copyright ©2024, Revvity, Inc. All rights reserved.